RANGER<sup>®</sup> 7000

**2<sup>nd</sup> Scale Interface**  
Instruction Manual

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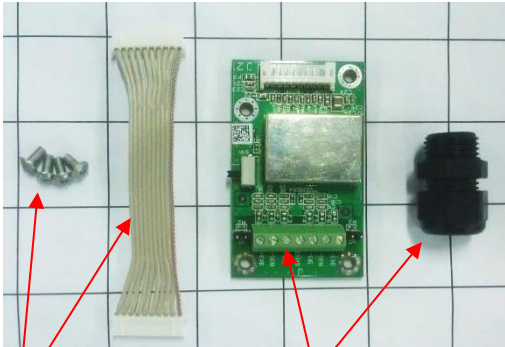
## INTRODUCTION

This Second Scale interface Kit is for use with the Ohaus Ranger<sup>®</sup> 7000.  
Through this kit you can connect an additional platform to the Ranger<sup>®</sup> 7000.

## KIT CONTENTS


- Screws (4)
- Cable to mainboard
- M16 Cable Gland
- 2<sup>nd</sup> Scale PC Board


## INTERFACE INSTALLATION

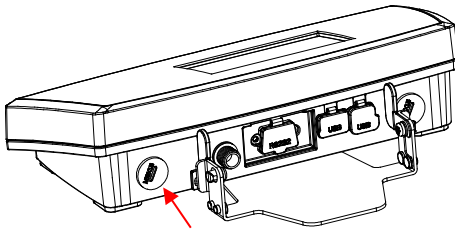Install the Interface module inside the terminal as shown.


1. Release the plastic plug by using a Slotted screwdriver

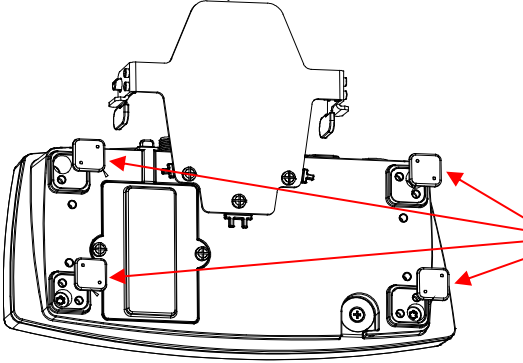
2. Remove the four rubber pads.

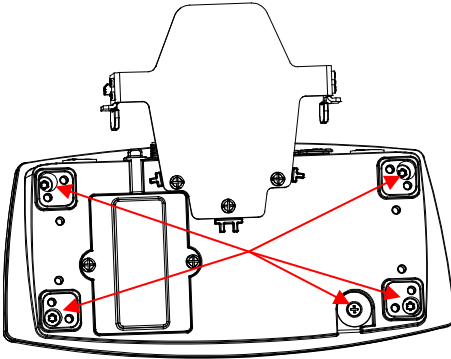
3. Release the five screws

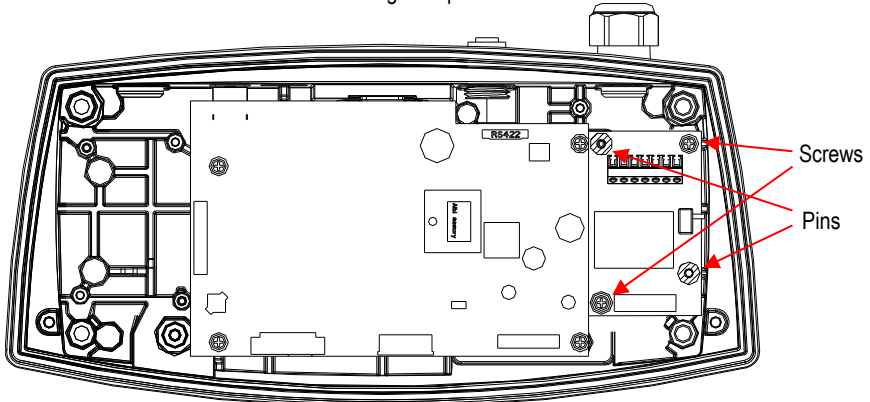
4. Install the Interface PCB. Push it through the pins and secure it with the 2 screws.

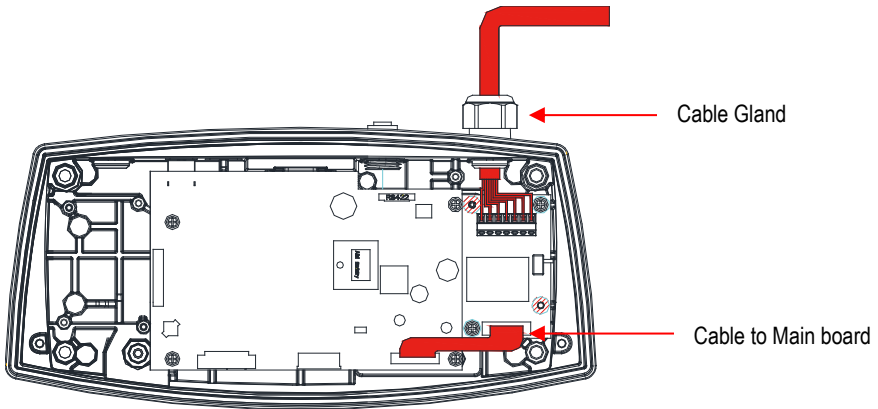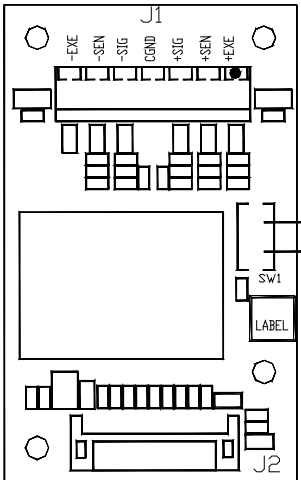5. Install the cable gland, and pass the load cell cable through it. Connect the cable between the A/D board and Main board.

6. Connect the load cell cable to the terminal block (J1).

7. Secure the cable gland and put the housing cover back on.

## SETUP

After connecting to the 2<sup>nd</sup> weighing platform, go to menu setup to set the 2<sup>nd</sup> Plat.Capacity, 2<sup>nd</sup> Plat.Cap.Unit and 2<sup>nd</sup> Plat.Grad and then perform a calibration.

Setup		
	Reset	>
	Auxiliary Display	On >
	Language	English >
	2nd Plat. Capacity	30
	2nd Plat. Cap. Unit	kilogram >
	2nd Plat. Grad.	0.001 >
		

### 2nd Plat.Capacity

Set the Capacity of Scale 2 using the numerical keypad

3 kg  
 ...  
**30 kg**  
 ...  
 10000 kg

### 2nd Plat.Cap.Unit

Use this menu to set the primary weighing unit for Scale2:

**Kilogram**  
 Pound

### 2nd Plat. Grad.

Use this menu to set the readability of scale 2

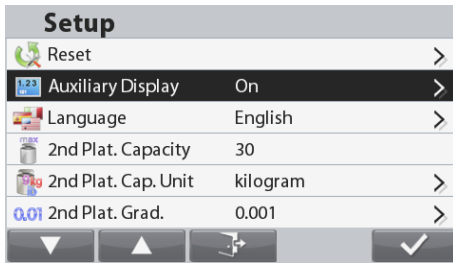
0.0005  
**0.001**  
 0.002  
 0.005  
 0.01  
 0.02

**Note:** When the scale2 is used in Legal for trade, Resolution <= 3000e.

Press and hold the  key on the overlay to switch between the main platform and the secondary platform.

**Note:** Make sure to perform a calibration of the scale before using the 2<sup>nd</sup> weighing platform.

To display the weighing results of both platforms, turn on the Auxiliary Display function in the SETUP menu (see menu structure below).

- Current platform icon
- Current platform weighing result
- Other platform weighing result

When the scale is used in trade or a legally controlled application, it must be set up, verified and sealed in accordance with local weights and measures regulations. It is the responsibility of the purchaser to ensure that all pertinent legal requirements are met.

Un-locked

Locked with Wire Seal

Locked with Paper Seal

**Note:** The Terminal only needs to be sealed if a second scale is attached to the optional 2<sup>nd</sup> A/D board.

## COMPLIANCE

### Disposal

In conformance with the European Directive 2002/96 EC on Waste Electrical and Electronic Equipment (WEEE) this device may not be disposed of in domestic waste. This also applies to countries outside the EU, per their specific requirements.

Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. If you have any questions, please contact the responsible authority or the distributor from which you purchased this device.

Should this device be passed on to other parties (for private or professional use), the content of this regulation must also be related.

For disposal instructions in Europe, refer to [www.ohaus.com/weee](http://www.ohaus.com/weee).

Thank you for your contribution to environmental protection.

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P/N 30105586 © 2013 Ohaus Corporation, all rights reserved / todos los derechos reservados / tous droits réservés / Alle Rechte vorbehalten / tutti i diritti riservati.
OHAUS Corporation