### Unpacking

Carefully unpack the Auxiliary Display kit. The package should contain:

- Auxiliary Display with attached cable
- Mounting brackets (2)
- This instruction sheet

### Installation

- Confirm the model number of your Auxiliary Display is appropriate for your balance (see table below).
- Unplug the balance.
- For TA, TAJ, SP, SPJ and SE Models: Connect the cable from the Auxiliary Display to the Option Port on the balance (Figures 1- 4). Re-connect power to the Balance.
- For PA, PAJ, PX and PR Models: Connect the RS232 and power cables from the Auxiliary Display to the receptacles at the rear of the balance (Figure 5). Then connect the power cable from the AC adapter to the receptacle on the cable of the Auxiliary Display. Re-connect power to the balance.
- Place the Auxiliary Display in the desired location.
- Turn on the balance and enter the balance's RS232 sub menu to configure the parameters as below:  
**Baud rate: 2400, Parity: 7 None, Stop Bit: 2, Handshake: None.**
- The Auxiliary Display operates automatically.

If communication is lost, turn the balance off and on to restore the connection with the Auxiliary Display.

### Display Mounting and Placement Options

Table top placement:

The Display may be placed on a flat surface without additional hardware.

Direct wall mounting:

The Display may be mounted to a wall using two screws (not included). Select appropriate sized screws that fit into the holes in the back of the Display.

Bracket wall mounting/table mounting:

Attach the two mounting brackets. Position the Display where desired on the wall or table.

Note: When installing the mounting brackets, insert round pins into holes in the back of the display and then slide each bracket to lock into place. Mounting brackets may be attached in two different orientations to provide optimum viewing angle.

### Technical Data

| Model                     | TAD7                                                                  | PAD7                                               |
|---------------------------|-----------------------------------------------------------------------|----------------------------------------------------|
| Ambient Temperature       | 5 °C to 40 °C                                                         |                                                    |
| Relative Humidity         | 15% to 80% at 31°C non-condensing, decreasing linearly to 50% at 40°C |                                                    |
| Height above Sea level    | Up to 2000m                                                           |                                                    |
| Storage temperature range | -10 °C to +60 °C                                                      |                                                    |
| Power requirements        | 3.3V === 30mA                                                         | 8-14.5V ~ 50/60Hz 40mA or 8-20V === 40mA (min-max) |
| Protection                | Pollution degree 2, Overvoltage category II<br>For indoor use only    |                                                    |

### Specifications

| Ohaus Balance         | TAD7 | PAD7 |
|-----------------------|------|------|
| TA, TAJ and SE Models | x    |      |
| PA, PAJ               |      | x    |
| SP, SPJ Models        | x    |      |
| EX, AX                |      | x    |
| PX, PJX               |      | x    |

This product complies with the applicable harmonized standards of EU Directives 2011/65/EU (RoHS) and 2014/30/EU (EMC). The EU Declaration of Conformity is available online at [www.ohaus.com/ce](http://www.ohaus.com/ce).

Register your accessory under *Warranty Registration at www.ohaus.com*

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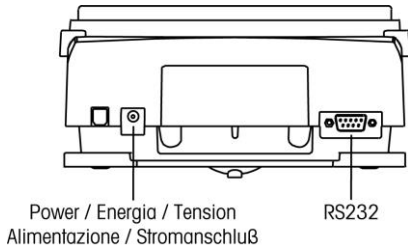
Figure 5 / Figura 5 / Figure 5 / Figura 5 / Abb 5

NOTE: This equipment has been tested and found to comply with the limits for a Class A digital device, pursuant to Part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment. This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this equipment in a residential area is likely to cause harmful interference in which case the user will be required to correct the interference at his own expense.

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LIMITED WARRANTY

Ohaus products are warranted against defects in materials and workmanship for duration of the warranty. During the warranty period Ohaus will repair, or, at its option, replace any component(s) that proves to be defective at no charge, provided that the product is returned, freight prepaid, to Ohaus. This warranty does not apply if the product has been damaged by accident or misuse, exposed to radioactive or corrosive materials, has foreign material penetrating to the inside of the product, or as a result of service or modification by other than Ohaus. In lieu of product registration, the warranty period shall begin on the date of shipment to the authorized dealer. No other express or implied warranty is given by Ohaus Corporation. Ohaus Corporation shall not be liable for any consequential damages. As warranty legislation differs from state to state and country to country, please contact Ohaus or your local Ohaus dealer for further details.

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P/N 80251398 D © 2016 Ohaus Corporation, all rights reserved / todos los derechos reservados / tous droits réservés / tutti i diritti riservati / Alle Rechte vorbehalten.